


## Assemble of the bioreactor




1: Put the blue silicon ring in the lid. Press down careful with a coin. 2: Put one silicon ring on the clamp.



3: Insert the clamp in one hole of the outer container and put the other silicon ring on the clamp from the other side and screw the nut on the clamp.

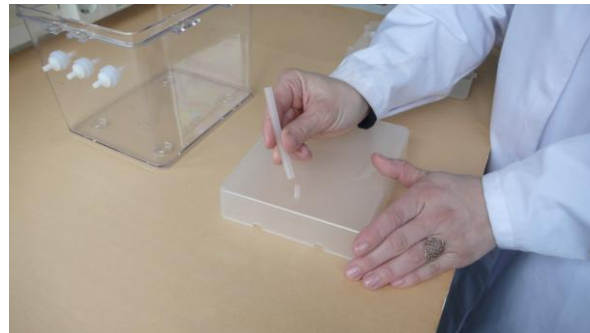

4: Do the same with the other clamps on the other holes. 5: Put the silicon tube on the inner chamber and put it in the bottom of the outer container.

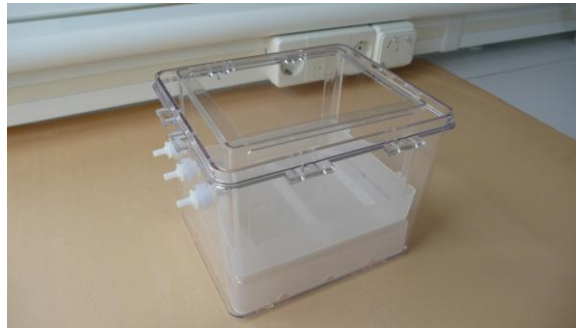

6: Put the basket above the inner container so the middle tube pass through the cut out of the basket.



7: Put the frame with the four legs above the basket and then close the container with the lid.